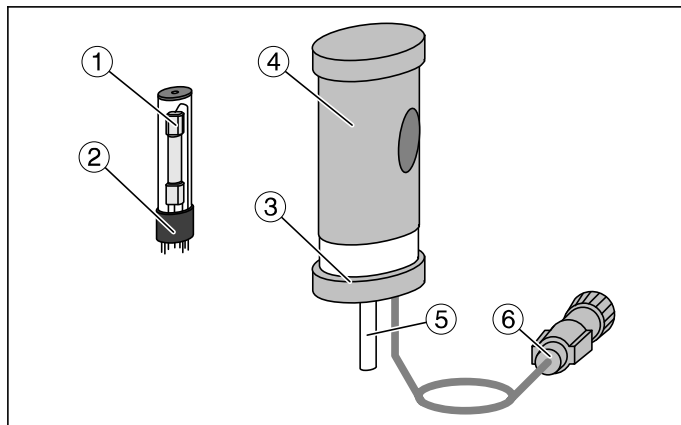
5/95-d.R.Sf-

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## Spectrum Lamps Ne, He, Cd, Hg, Hg-Cd, Tl, Na Housing for Spectrum Lamps

Fig. 1

The spectrum lamps are AC-operated gas-discharge lamps emitting the line spectrum of metal vapors or noble gases respectively with high spectral purity and high luminous intensity.

The spectrum lamps are set up in the housing for spectrum lamps (451 16); the power lead of the housing is connected to the universal choke (451 30), which supplies precisely the current needed to operate the lamps.

#### Examples of Experiments:

- Determination of wavelengths
- Inversion of the Na-D line (using 451 111)
- Investigation of UV radiation (using 451 062)
- Photoelectric effect (using 451 062)
- Determination of  $h$  (using 451 062)

## 1 Safety notes

Spectral lamps (451 011 and following numbers) in housing (451 16) should be connected to the mains only via the universal choke (451 30)

Keep a minimum distance of 3 cm between light exit opening and optical element (e.g. diaphragm, lens) to avoid overheating due to accumulation of heat.

## 2 Technical Data

### 2.1 Spectrum lamps (451 011 to 111)

- ① Glass bulb (10 cm x 2 cm dia.) with burner
- ② 9-pin pico socket matched to housing (451 16)

Gas filling*)	Ne	He	Cd	Hg	Hg-Cd	Tl	Na
Cat. No.	451 011	451 031	451 041	451 062	451 071	451 081	451 111
Operating voltage	30 V	60 V	15 V	45 V	30 V	15 V	15 V

Operating current:  $1 \text{ A} \pm 10\%$

(ensured with supply via universal choke, 451 30, at mains voltages between 200 V AC and 230 V AC)

\*) Spectrum lamps filled with caesium, potassium, rubidium and zinc are also available to special order.

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## 2.2 Housing for spectrum lamps (451 16)

- ③ Lower part of housing with 9-pin pico socket, built-in starter
- ④ Upper part of housing, movable on ③, with light emission aperture (2.5 cm dia.)  
Dimensions of housing: 16 cm x 7 cm dia.
- ⑤ Stand rod, 5.5 cm long, 10 cm dia.
- ⑥ Connecting lead, 80 cm long, with multipin plug, to be connected, safe to touch, to the universal choke (451 30) by means of union nut.

## 3 Operation

*Additionally required:*

Universal choke 451 30  
Optical elements and  
Stand material, e.g. according to Fig. 2

Plug in spectrum lamp, connect housing with universal choke, switch on.

Allow to warm up, the full radiation output is achieved only after some minutes, except spectrum lamp He (451 031) and spectrum lamp Ne (451 011).

In case of power failure, allow the warm spectrum lamp Hg (451 062) to cool down some minutes before switching it on again (otherwise it will not ignite).

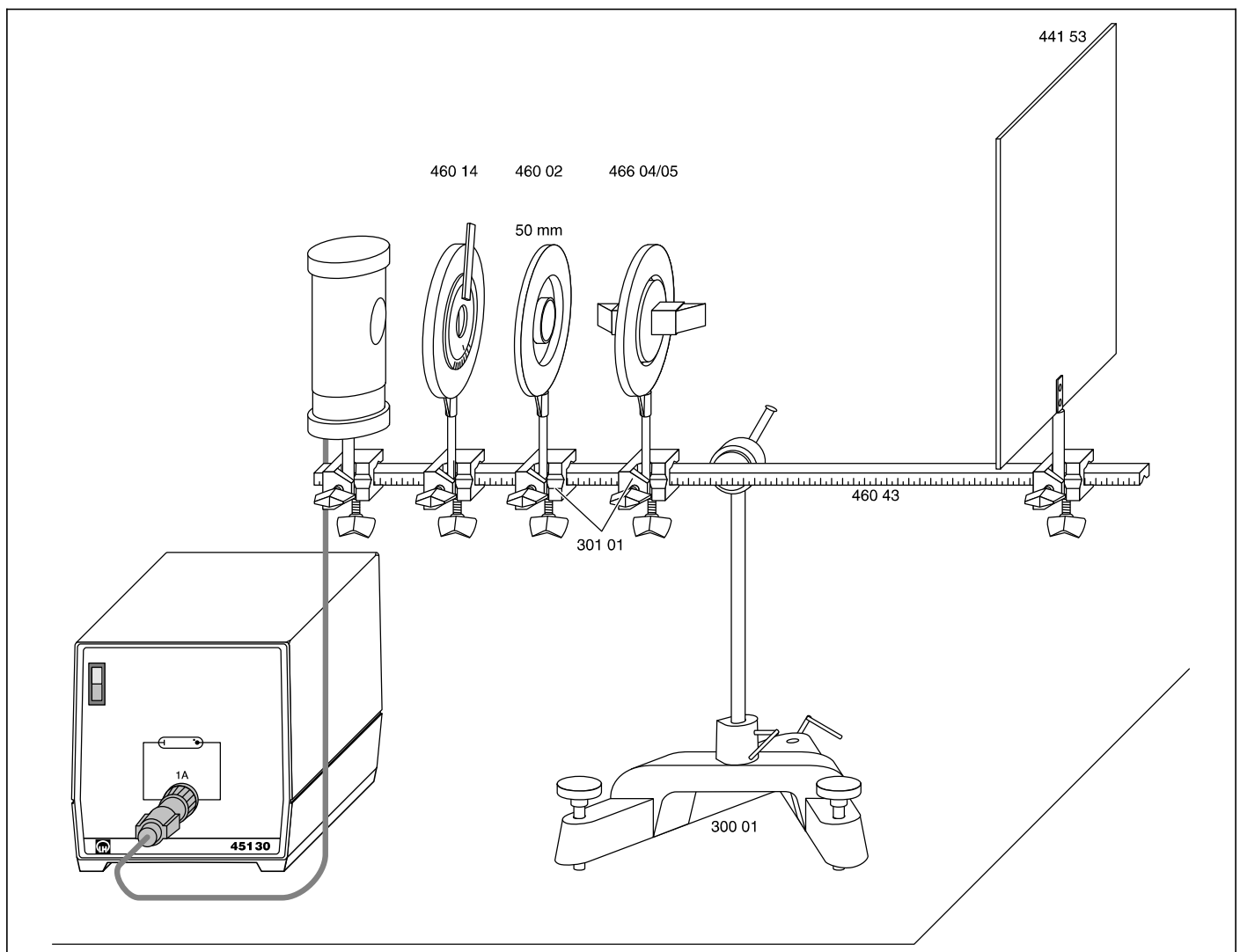
Fig. 2