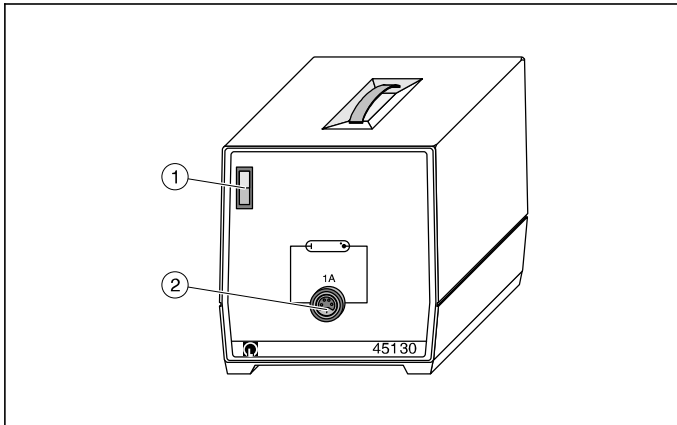
3/96-Sf-

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451 30*

Universaldrossel in Gehäuse Universal Choke in Housing

Fig. 1

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The universal choke is designed for use with 230 V/50 Hz voltage for supplying (1 A) the apparatus:

- High-pressure mercury lamp (451 15) in lamp socket E 27 (451 19, series 3 on) or in the compact arrangement for determining Planck's constant (558 79, series 2 on)
- Cadmium lamp for Zeemann effect (451 12), and
- Spectral lamps (451 011 - 451 111) in housing (451 16, series 2 on).

1 Safety notes

- Do not switch on the apparatus until the special plug of the socket in which the lamp is mounted has been correctly connected to the DIN socket ② (secure the connection using the union nut of the special plug).
- Apparatus from older series (than those specified above) must be returned to Leybold Didactic for conversion to match the universal choke (451 30, series 4).

2 Description, technical data

- ① Mains switch with integrated indicator lamp
- ② DIN socket
Output current: 1 A

On rear of housing:

Appliance plug connector for mains power lead (included in scope of supply)

Fuse holder, integrated in appliance plug, with operating and spare fuse T 1.25 B

Mains connection voltage: 230 V/50 Hz
Dimensions: 20 cm x 21 cm x 23 cm
Weight: 5 kg

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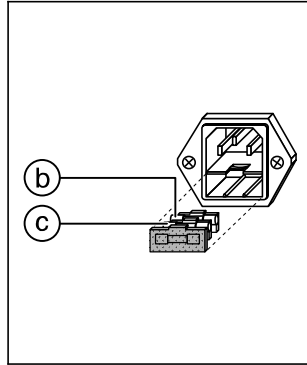
Fig. 2.2

3 Replacing the fuse (see Fig. 2.1/2)

- Pry out insert (a) (with holder for primary fuse (b) and reserve fuse (c)).
- Replace defective fuse (b) with a new one which has been checked for the correct rating (c).
- Insert reserve fuse (c) and replace the insert (a) (Cat. No. for 10 fuses T 1,25 B: 698 16).