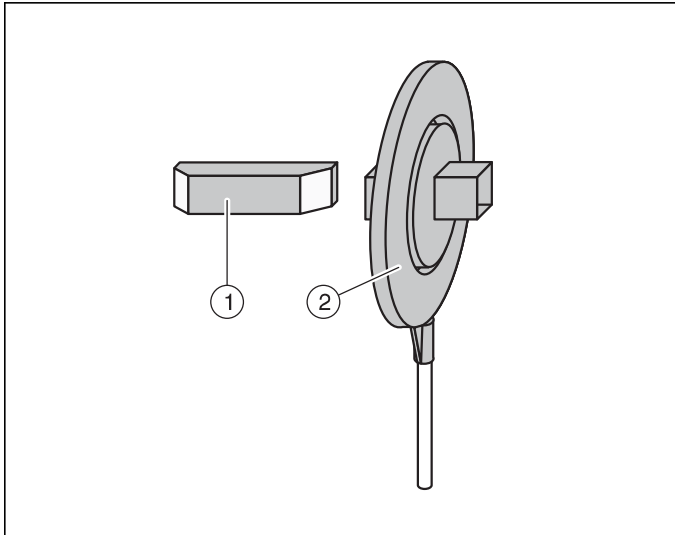
9/97-Sf-

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The direct vision prism allows spectral dispersion of light without changing the direction of the beam axis.

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Examples of experiments

- Dispersion and recombination of white light; complementary colours
- Absorption spectra (coloured glasses, liquids, gases; band and line spectra)
- Emission spectra (line spectra of inert gases and metal vapours)
- Reversal of the Na-D-line
- Fraunhofer lines
- h -determination from the Hg spectrum

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1 Safety notes

When inserting the prism into the holder, do not toe it (risk of breakage).

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2 Description, Technical Data

- ① Direct vision prism (466 05),
made up of 3 prisms (1 x flint glass, 2 x crown glass)
Angular dispersion (C - F): $4^{\circ} 14'$
Cross-section: 20 mm x 20 mm
Length: 10 cm
- ② Holder for direct vision prism (466 04)
Prism table, eccentrically arranged so that when fitting the holder to an optical bench the refraction surface of the prism coincides with the optical axis.
On stand rod (10 cm length, 10 mm dia.).

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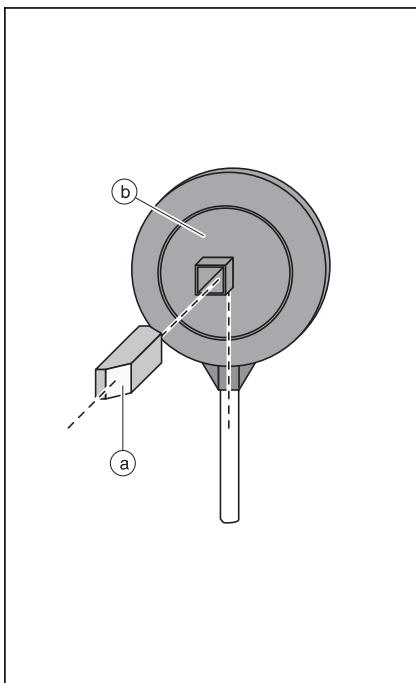
Fig. 2

3 Use

Prior to first-time operation, push the prism into the holder as shown in Fig. 2 so that the refraction surface (a) is in the middle of the holder surface (b).

Important: Do not toe the prism when inserting it into the holder (risk of breakage).

If due to manufacturing tolerances the prism should not fit into the holder without slipping, secure it by adhesive tape.

Additionally required for optical assembly (see Fig. 3)

Source of light according to experimental conditions, e.g.

Lamp housing with lamp, 6 V, 30 W	450 60/51
Lamp housing with halogen lamp, 12 V/100 W	450 63/64
Spectrum lamps with housing	451 011-111/16
High-pressure mercury lamp with socket E 27	451 15/19

Condenser, e.g.

Aspherical condenser for 450 60	from 460 20
Lens in holder, $f = 50$ mm	460 02

Adjustable slit 460 14

Imaging lenses, e.g.

Lens in holder, $f = 100$ mm	460 03
or	
Projection objective, $f = 150$ mm	460 13

Translucent screen 441 53

Optical bench with riders, e.g.

Small optical bench with Leybold multiclamps	460 43/301 01
---	---------------

Focus the light through the condenser onto the slit and form a sharp image of the slit on the screen before placing the direct vision prism into the beam of light.

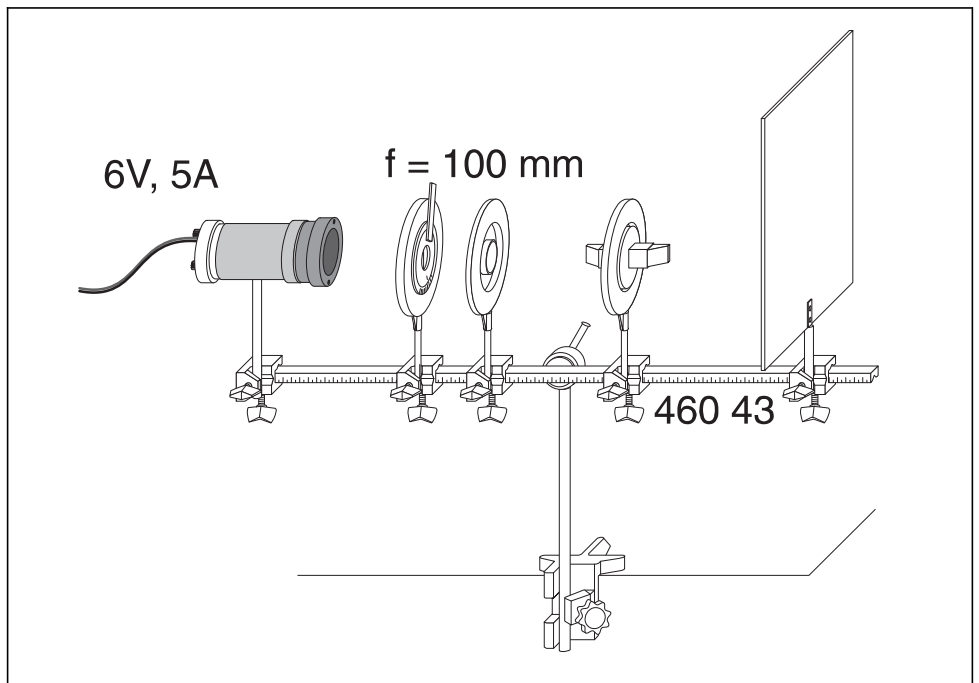
Fig. 3